



Catalog Number:	C1010
Cell Line Name:	CHO suspension cell pool expressing full-length human LAG-3 receptor
Gene name:	LAG-3, human
Gene Sequence:	Codon optimized from NP_002277.4 (Met1-Leu525)
Host Cell:	Suspension CHO
Quantity:	Two vials of frozen cells (2×10^7 per vial)
Lot Number:	12042
Stability:	>10 passages
Freeze Media:	Culture media with 10% DMSO
Storage:	liquid nitrogen immediately upon receipt
Culture Medium:	50% CD-CHO (Gibco [#] 10743-029), 50% Ex-Cell CHO 5 Media (Sigma [#] C0363), supplemented with 8mM L-Glutamine, 1xHT, 1x Penn-Strep and 20ug/ml puromycin.
Mycoplasma Test:	Negative
Application:	Antibody binding assays, IHC/Western blot analysis, or use as cell immunogen

I. Background:


Lymphocyte-activation protein 3 (LAG-3) belongs to the immunoglobulin superfamily and shares sequence homology, exon/intron organization and chromosomal localization to CD4. LAG-3 cDNA encodes a 498 amino acid transmembrane protein with 4 extracellular Ig-like domains. The expression of LAG-3 is undetectable in resting peripheral blood lymphocytes but it is induced in activated T and NK cells. Human MHC Class II molecules can bind to LAG-3 and are considered as receptors for LAG-3 on B cells and dendritic cells. LAG-3 is proven to be an inhibitory receptor on activated T cells. Crosslinking of LAG-3 on activated T cells with an anti-LAG-3 antibody inhibits T cell proliferation and cytokine secretion. Anti-LAG-3 antibodies which blocked LAG-3 binding to MHC class II molecules had strong anti-tumor activity when used in combination with anti-PD-1 antibodies in tumor xerograph models.

Disclaimer: For research use only. Not for use in humans.



II. Representative Data:

Detection of human LAG-3 expression on human LAG-3 CHO Stable cells using a mouse monoclonal antibody specific for human LAG-3.

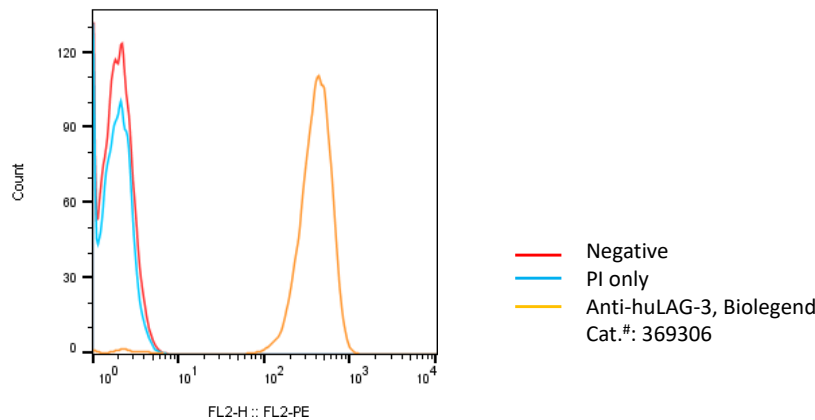

III. Thawing and Subculturing:

1. Remove the cell vial from liquid nitrogen tank and thaw cells quickly in a 37°C water bath.
2. In a T250 shake flask, add 30ml cell culture media which has been pre-warmed in a 37°C water bath for 30 minutes.
3. Transfer all the cells in the vial into the T250 shake flask and place it on a shaker inside of a CO₂ incubator with 5-8% CO₂. The shaker speed should be set at 120-150rpm.
4. The next day, count the cells using a hemocytometer. The viable cell density should be around 0.5-1x10⁶/ml and the cell viability should be more than 50%.
5. The cell doubling time should be 18-24 hours. Split the cells 1:5 or 1:10 every 2-3 days or when the cell density reached 6-8x10⁶/ml.

IV. References:

Triebel F. et al., LAG-3, a novel lymphocyte activation gene closely related to CD4. *J Exp Med*, 1990 May 1; 171(5): 1393-405.

Baixeras E. et al., Characterization of the lymphocyte activation gene 3-encoded protein. A new ligand for human leukocyte antigen class II antigens. *J Exp Med*, 1992 Aug 1;176(2):327-37.



Hannier E., et al., CD3/TCR complex-associated lymphocyte activation gene-3 molecules inhibit CD3/TCR signaling. J Immunol. 1998 Oct 15;161(8):4058-65.

V. Limited Usage License Agreement:

This is a legal agreement between the buyer (Licensee) and Accurus Biosciences Inc. regarding the use of Accurus Biosciences' stable cell line products. By purchasing and using the stable cell lines, the buyer agrees to comply with the following terms and conditions of this label license and recognizes and agrees to such restrictions:

- A) The products are not transferable and will be used at the site where they were purchased. Transfer to another site owner by the buyer or to a subcontractor will be permitted only when written approval from Accurus Biosciences is obtained.
- B) The buyer cannot sell or otherwise transfer a) this product, b) its components or c) materials made using this products to a third party.
- C) The products sold by Accurus Biosciences are for laboratory and animal research use only. The products are not to be used on human studies or consumption.

Disclaimer: For research use only. Not for use in humans.